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Register for a JEOL Delta/NMR Support Site Account

http://nmrsupport.jeol.com/Home/Register/tabid/53/Default.asp

LogIn Register

Wednesday, November 10, 2010

**JEOL**  
NMR SUPPORT SITE

Home Go

### JEOL Delta/NMR Support Site Account Registration

Completely fill out this form. Your request will be forwarded to the site administrator for review and acceptance. You will be contacted once your account has been approved.

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
The screenshot shows a web browser window titled "Register Spectrometer" with the URL "http://nmrsupport.jeol.com/Home/RegisterSpectrometer/tabid/77". The page header includes the JEOL logo and "NMR SUPPORT SITE". The main content area is titled "Spectrometer and System Information" and contains a form with the following fields and annotations:

- Spectrometer Brand:** JEOL (dropdown menu) - Annotation: • NMR装置メーカー
- JEOL Spectrometer Type:** ECA (dropdown menu) - Annotation: • NMRの機種
- Spectrometer Field:** 600 (dropdown menu) - Annotation: • 磁場強度
- Workstation:** Intel PC (dropdown menu) - Annotation: • PCの種類
- Operating System:** Windows (dropdown menu) - Annotation: • OSの種類
- Operating System Version:** XP Professional (dropdown menu) - Annotation: • OSのバージョン
- Do you need Delta for:** Radio buttons for "Data Processing Only" (selected), "Data Processing and JEOL NMR Machine Control (JEOL NMR Spectrometers Only)", and "None". Annotation: • 処理用を選択します
- Are you a JEOL employee or representative?:** Radio buttons for "No" (selected) and "Yes". Annotation: • Noを選択します
- Submit:** A button to submit the form. Annotation: • 全て入力したらクリックします

Additional text on the page includes: "Your registration is **NOT** complete until you supply the information requested below.", "Logout Katsuo Asakura", "Thursday, November 11, 2010", and contact information for JEOL USA, Inc. at the bottom.

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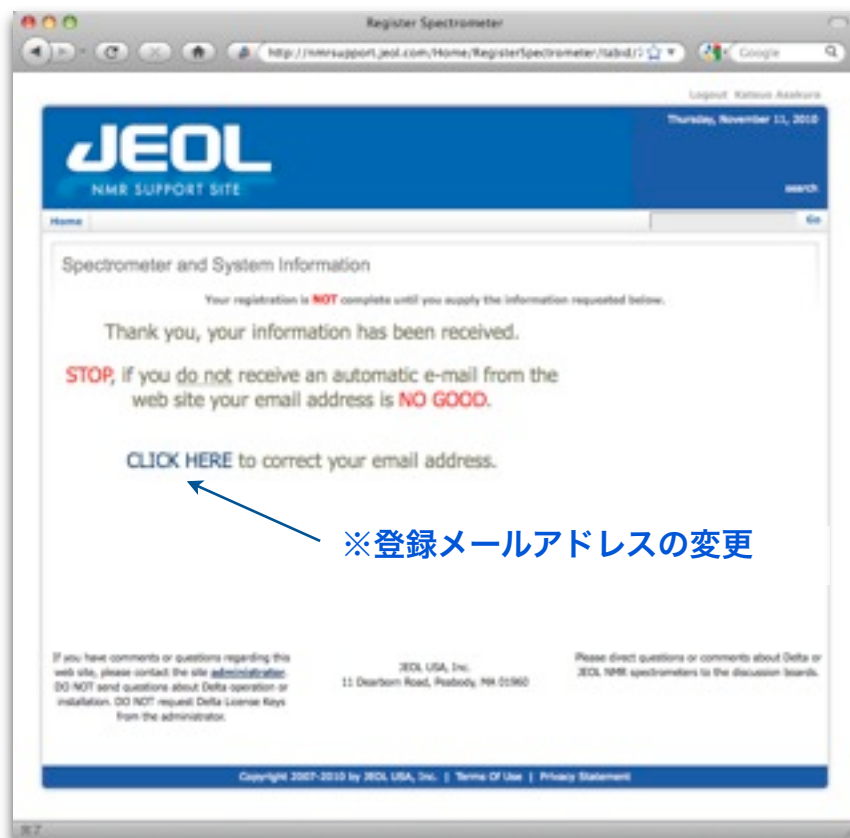Logout Katsuo Asakura  
Thursday, November 11, 2010

**JEOL**  
NMR SUPPORT SITE

Home Software Calendar Docs Forums FAQs Go

Welcome to JEOL Delta / NMR Support

The JEOL Delta / NMR support web site provides support for users interested in the Delta NMR Software package and JEOL's NMR Spectrometers. The web site includes Delta NMR software kits for NMR data processing only and kits for JEOL NMR Spectrometer control with data processing. Areas for Delta License Key generation, discussion forums, FAQ's, and documentation downloads are also available on this web site.

To get started, simply:

- Download Delta Software** ← ダウンロードページへ  
Depending on your account type, Delta is available for data processing only, or for full JEOL NMR spectrometer control. The current and older versions of Delta are available for download.
- Obtain a License Key**  
Submit the Delta Machine ID of your workstation to obtain a matching Delta License Key to enable the software.
- Document Downloads**  
Download Release Notes, Delta Tutorials, User Guides, and Other Documents
- Discussion Forum**  
Discussion Forums Related to Delta and JEOL NMR Spectrometers
- FAQs**  
Frequently Asked Questions related to Delta and JEOL NMR Spectrometers.
- Manage Profile / Change Password**  
Manage contact and spectrometer information or change password. The spectrometer and workstation information are on the second page of the profile after "UPDATE" of the first page.

Read Me 日本語

Title  
NMR Support Read Me 日本語 (Japanese)  
NMR Support Read Me (English)

Get ADOBE READER

If you have comments or questions regarding this web site, please contact the site [administrator](#). DO NOT send questions about Delta operation or installation. DO NOT request Delta License Keys from the administrator.

JEOL USA, Inc.  
11 Dearborn Road, Peabody, MA 01960

Please direct questions or comments about Delta or JEOL NMR spectrometers to the discussion boards.

Copyright 2007-2010 by JEOL USA, Inc. | [Terms Of Use](#) | [Privacy Statement](#)

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Software

http://nmrsupport.jeol.com/Software/tabid/75/Default.aspx

Logout: Katsuo Asakura  
Thursday, November 11, 2010

**JEOL**  
NMR SUPPORT SITE

Home Software Calendar Docs Forums FAQs

### Delta Software Download

#### Terms and Conditions

CAREFULLY READ THE FOLLOWING TERMS AND CONDITIONS. YOU AGREE TO ALL TERMS AND CONDITIONS BY INSTALLING THE DELTA NMR SOFTWARE PACKAGE OR USING ANY PART THERE OF.

License: With the payment of any and all applicable License Fees upon the delivery of Delta NMR Software ('Software'), JEOL USA, INC. ('Licensor') grants to the Licensee a License to use without exclusivity a single non-transferable, non-sublicensable copy of the

Accept Terms  Reject Terms

Package Type:  Data Processing Only  Full Package with Instrument Control

Workstation: Intel PC

Operating System: Windows

Operating System Version: XP Professional

Files to Review & Retrieve:  Current Version Only  All Files

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Available Software Downloads (all versions)

Delta for Windows / v4.3.6

Show Available Software Download Software

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License Keys

http://nmrsupport.jeol.com/LicenseKeys/tabid/76/Default.aspx

Logout Katsuo Asakura  
Sunday, November 21, 2010

**JEOL**  
NMR SUPPORT SITE

Home Software License Keys Docs Forums FAQs

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Accept Terms  Reject Terms

Data Processing Only  Full Key with Instrument Control

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If you have comments or questions regarding this web site, please contact the site administrator. DO NOT send questions about Delta operation or installation. DO NOT request Delta License Keys from the administrator.

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Copyright 2007-2010 by JEOL USA, Inc. | Terms Of Use | Privacy Statement

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License Keys

http://nmrsupport.jeol.com/LicenseKeys/tabid/76/Default.aspx

Logout Katsuo Asakura  
Sunday, November 21, 2010

# JEOL

NMR SUPPORT SITE

Home Software License Keys Docs Forums FAQs

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Accept Terms  Reject Terms

Data Processing Only  
 Full Key with Instrument Control

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Delta Machine ID: SJ - VH - SO - WN - HX

Your Delta Machine ID is SJ-VH-SO-WN-XXXX

Your Delta License Key is HS-TS-XV-PQ-XXXX

This information is also being E-mailed to your address on record.

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If you have comments or questions regarding this web site, please contact the site administrator. DO NOT send questions about Delta operation or installation. DO NOT request Delta License Keys.

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